

Product Datasheet

STD-68G Water Treatment Additive



An 8 fl oz (236ml) bottle of water preservative for use with the STD-68G portable self-contained eye/face wash stations. It preserves from 15-76 litres of water and two bottles come free when purchasing an STD-68G eye/face wash station. The station should be cleaned and refilled every 120 days when using the water preservative.

Specifications

INGREDIENTS

Chlorhexidine 1% w/v Purified Water

Propylene Glycol

USAGE

For use with the STD-68G self-contained emergency eye wash station

WARNINGS

For external use only. In case of contact with eye in undiluted form, flush with clean water

WHEN USING THIS PRODUCT

Do not change dilution or use with other chemicals

Do not reuse after application

SEEK MEDICAL ADVICE IF YOU HAVE ANY OF THE FOLLOWING DURING OR AFTER USE:

Changes in vision

Eye pain

Condition worsens or persists

Continued redness or irritation of the eye

If swallowed

DIRECTIONS

Wear protective eyewear and gloves

Clean STD-68G and rinse with potable water

Partially fill STD-68G with potable water

Remove tamper evident seal and cap of bottle

Add entire contents of the bottle to the STD-68G

Fill the STD-68G to the required fill level

Date and initial inspection tag

STD-68G should be cleaned and refilled every 120 days

OTHER INFORMATION

Batch number is printed on the bottle

Store at 20°C - 25°C

Do not use if seal is missing or broken

Use before expiration date marked on the bottle

Once opened, use entire contents and discard

Expiration date only applied to a sealed bottle

Keep out of reach of children

View our Guide to the ANSI/ISEA Z358.1-2014 International Standard for Emergency Eyewash and Shower Equipment at www.hughes-safety.com



HUGHES



HUGHES-SAFETY.COM

Justrite Safety Group

www.hughes-safety.com